

## **UD Cohesive Sampler**

## Unit Dose Sampler for Cohesive Powders

A popular sampler in the pharmaceutical industry for taking low volume (unit dose) cohesive powder samples. Allows accurate samples to be taken.

Type: Single Point Unit Dose

Material: 316 stainless steel

Operation: 1. Insert the sampler into the product.

- At the required depth push down the handle and rotate it through 180°.
- Pull up sampler several millimetres (to fill the cell), then rotate handle 180° (to lock cell in closed position).
- 4. Remove sampler




A) 'T' Bar Handle

Special lengths available on request

C) Polished finish

Sample
Collecting Cell different sized
cells can be
easily
interchanged

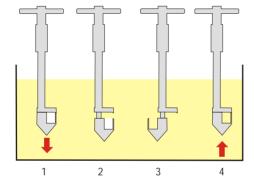



'T'Bar Handle



Tip in open position



Tip in closed position









Part No.	Description
2000S-01	U-D Cohesive Sampler Set - consisting of:
	1 off Handle, Sample Collecting Cells - 1 each of: 0.25ml, 0.5ml, 0.75ml and 1.0ml,
	1 off Superior Sampler Case
2000A-1000	Sampler body and handle - 1000mm length
2000T-025	Sample Collecting Cell - 0.25ml
2000T-050	Sample Collecting Cell - 0.5ml
2000T-075	Sample Collecting Cell - 0.75ml
2000T-100	Sample Collecting Cell - 1.0ml
8001N-1100	Cleaning Brush - 1100mm length - narrow style
	Storage Cases - download Range Brochure for further information