

## **U-D Master**

## Unit Dose Sampling to meet FDA Guidelines


Designed for multilevel low volume sampling in the Pharmaceutical industry. It is ideal for use with free flowing powders. By simply changing the sample inserts (cells) different volumes can be sampled.

More than one sample can be taken from the same sampling point (see double or triple styles).

Material: 316 stainless steel, satinfinish


Operation:

- Select the correct volume sampling cells and fit them to the inner core.
- 2. Locate the assembled inner core into the outer sleeve at the lowest sampling level
- 3. Insert the sampler into the product, turn the handle to expose the sampling cells to the product.
- Double and Triple Style samplers pull the handle to the next level to expose the next set of sampling cells.
- Turn the handle so no sampling inserts are exposed then remove the sampler from the product.
- Remove the inner core from the outer sleeve, the individual sampling cells can then be easily removed for analysis.

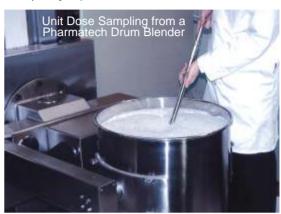




Place the Sampling cells in the inner core

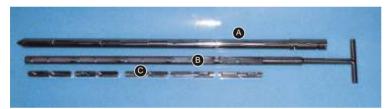

- (A) Outer Sleeve
- (B) Inner Core

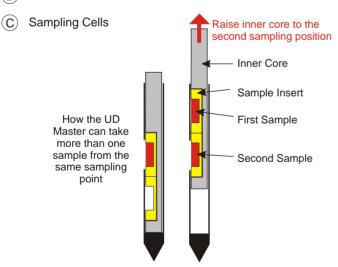

SINGLE Style Sampler

Collects ONE set of samples at each sampling point.

**DOUBLE** Style Sampler

Collects TWO sets of samples at each sampling point.

TRIPLE Style Sampler

Collects THREE sets of samples at each sampling point.



Sampling Cells are stamped with their volume for easy identification



Bayonet fitting allows two samples to be taken at the same sampling point



Slide the inner core into the outer sleeve

Tel: +44 (0)1675 466992

Fax: +44 (0)1675 466994

www.samplingsystems.com

e-mail: info@samplingsystems.com



## **U-D Master cont....**









	Part No.	Sampler Length	Number of Sampling Points	Number of Sample Cells needed
	1230A-300	300mm	5	5
SINGLE	1230A-600	600mm	10	10
	1230A-900	900mm	6	6
Style Sampler	1230A-1200	1200mm	8	8
	1230A-1500	1500mm	10	10
	1230A-1800	1800mm	12	12
DOLIDI E	1230B-600	600mm	5	10
DOUBLE	1230B-900	900mm	7	14
Style Sampler	1230B-1200	1200mm	8	16
21,712 23.11.12.12.1	1230B-1500	1500mm	10	20
	1230B-1800	1800mm	12	24
	_			
TRIPLE	1230C-900	900mm	5	15
Style Sampler	1230C-1200	1200mm	7	21
Style Sampler	1230C-1500	1500mm	8	24
	1230C-1800	1800mm	10	30
Sample Cells	S -			
Part Number is 12:	vailable in 0.1ml increments 30T-XX (where XX = Cell V	olume)		
For Example - The Part Numbers for 1.2ml and 0.5ml cells are: 1230T-12 and 1230T-05 respectively				
Accessories	Cleaning Brushes	Download Range Brochure for full details		
	Storage Case	Download Range Brochure for full details		